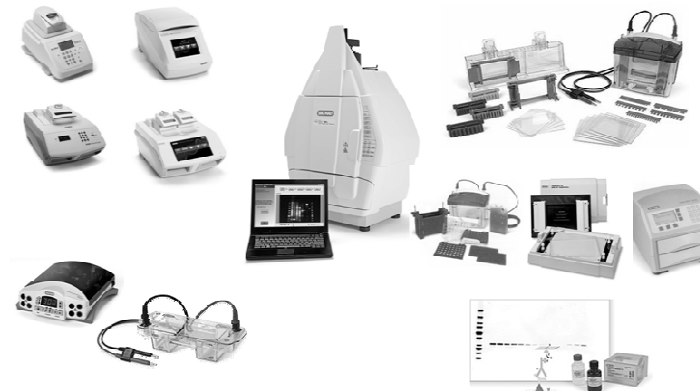
Imaging System Geldoc XR⁺

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
Applications Testing

- Applications specific testing for camera validation:
- DNA Applications:
 - EtBR = 100 pg sensitivity
 - SYBR Green = 50 pg
- Protein Applications:
 - Oriole < 1ng
- Bio-Safe Coomassie = 4 ng

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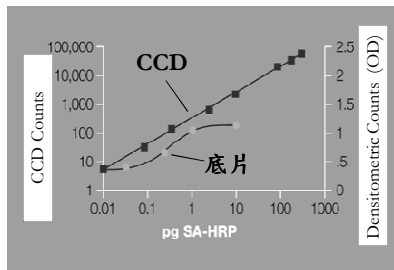
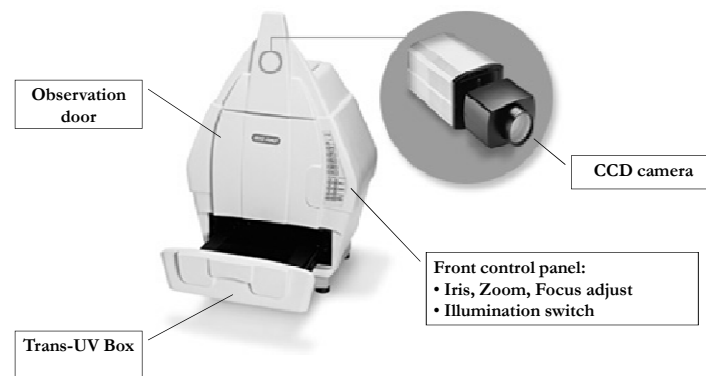
GelDoc XR+

GelDoc XRS+
-Configurations

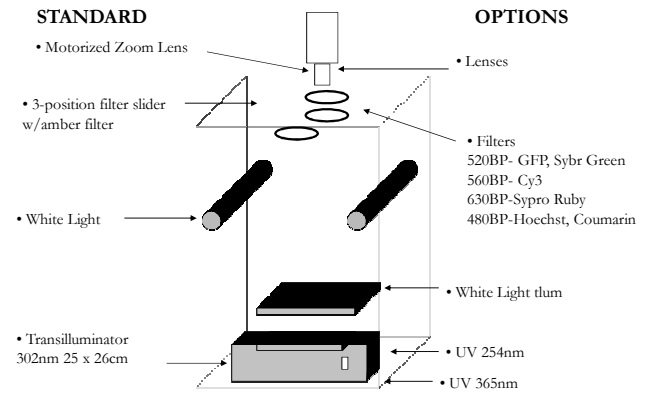
Unique Features – Dynamic Flat Fielding

- Patent-pending Bio-Rad technology
- Elimination non-uniformity caused by variations in the optical path due to zoom setting, illumination, and/or lens.
- Demonstrates the excellent uniformity of the system

Tech Note #2753

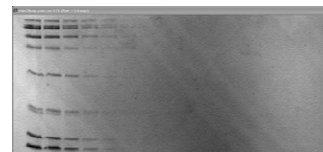Benefit of Flat Fielding

No flat field

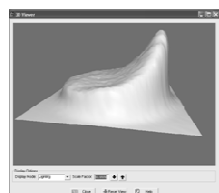
With flat field

BioSafe Coomassie stain was incubated overnight on a 8-16% Tris-HCl gel to detect Broad Range markers. Flat fielding removes background non-uniformity.

VersaDoc MP

GelDoc XR Comparison


Slightly saturated Harta LED spot – displayed as single peak

Spot #1 – Slightly Saturated

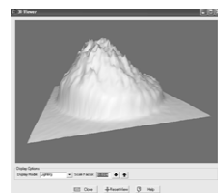
Un-saturated Harta LED Spot – displayed with rounded peak

Spot #3 – Not Saturated

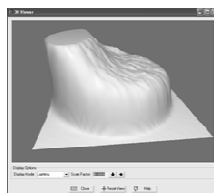Over-saturated Harta LED Spot – Displayed with flat peak

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Introduction to Software Image Lab

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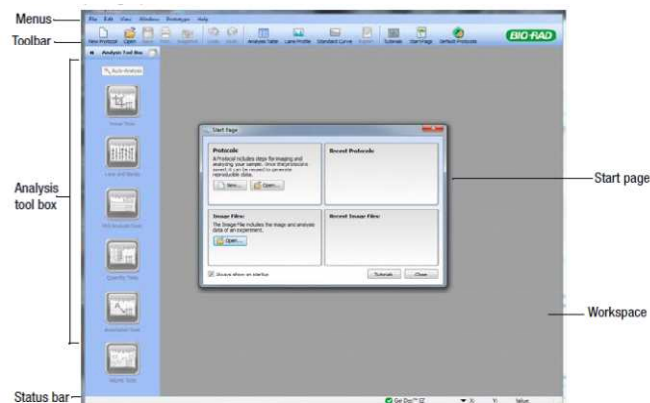
Easy Software Include

- **FAST**- One step imaging and analysis
- **Automatic Iris, Focus**
- **Resolution**
- **Reproducibility**
- **Saturation display**
- **Report**

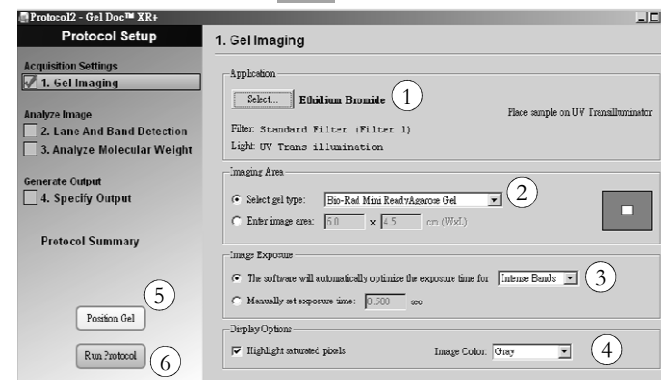
Image Lab Software

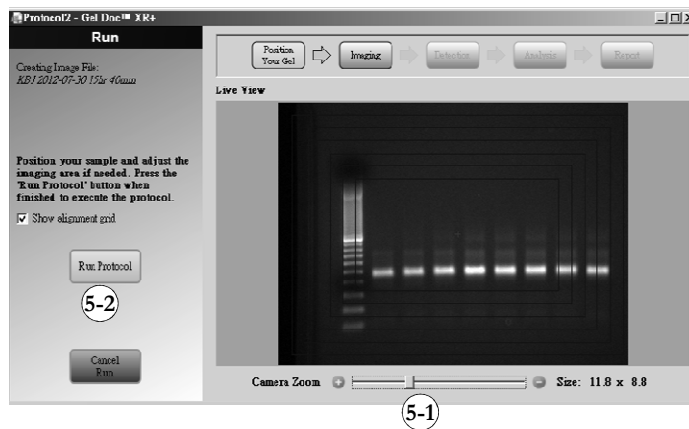
Protocol for imaging blot on GelDoc XRS+ using Image Lab software


- Click on the new protocol icon

Step 5, Check gel size and zoom in/out

Get result and save it.

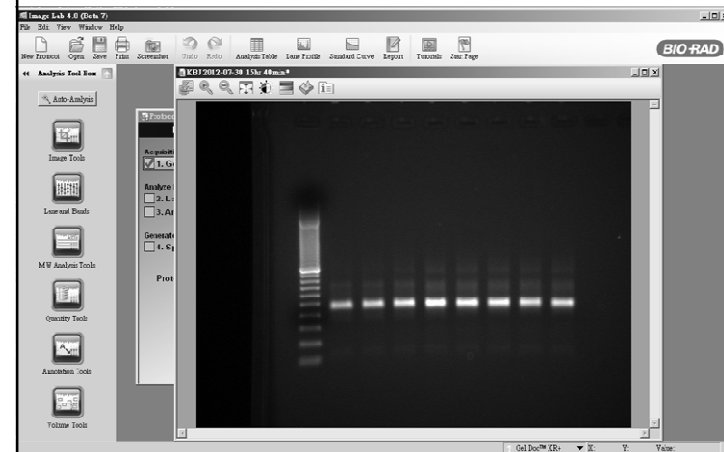Export TIFF or Jpeg

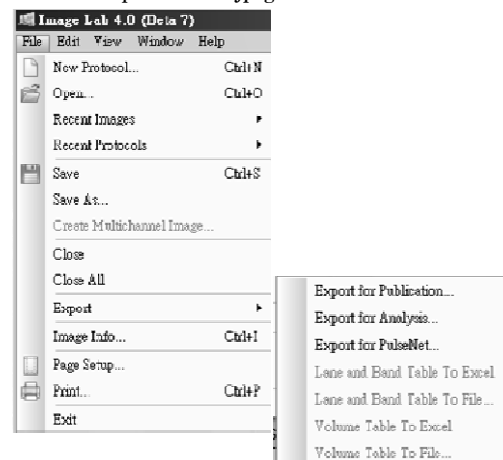
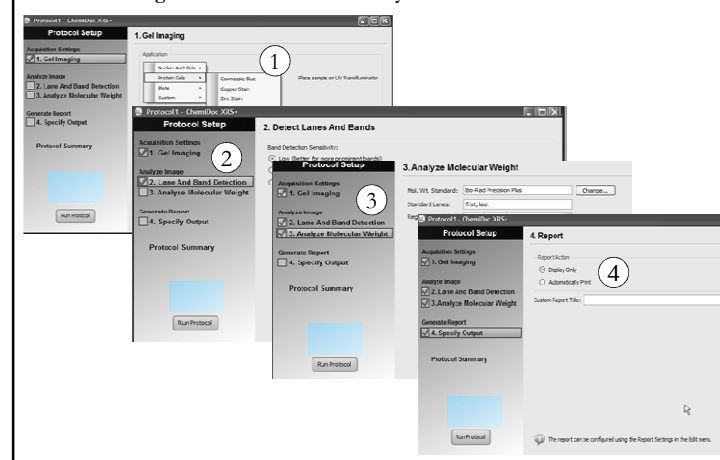
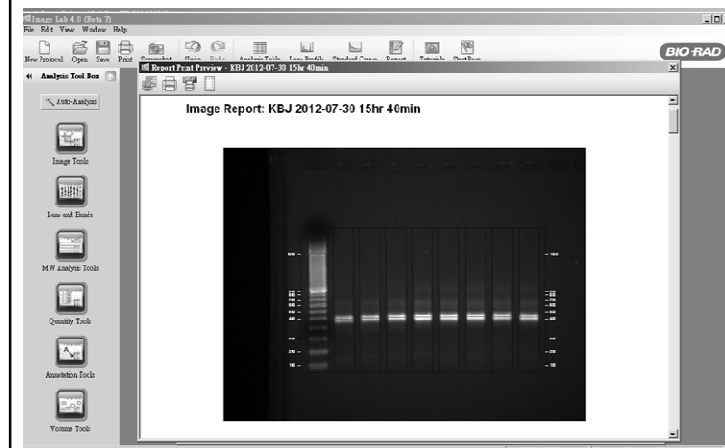
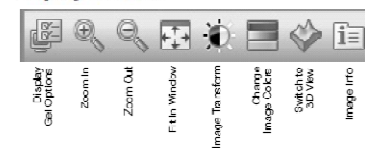
Image Lab – Get result and analysis it in the same time

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